

ELISA Accessory Kit

Cat#:1030011

[Product Description]

ELISA auxiliary reagent contains 8 main auxiliary components in ELISA kit, excluding antibody, enzyme, standard and other control substances. Please check the composition of the reagent before use. If you have any questions, please contact Shenzhen Dakewe Bio-engineering Co., Ltd.

[Product Composition]

Product Name	Cat No.	Component Name	Size	Quantity	Reconstitution
ELISA Auxiliary Kit Manual	1030011	Coating Buffer	12 mL	1 bottle	Ready-to-use, 100 μ L/well
		Blocking Diluent	25 mL	1 bottle	Ready-to-use, 250 μ L/well
		TMB	10 mL	1 bottle	Ready-to-use, 100 μ L/well
		Dilution Buffer R(1 \times)	10 mL	3 bottles	Ready-to-use
		Washing Buffer (50 \times)	15 mL	1 bottle	Dilute 1:50 with distilled water into working solution, 300 μ L/well
		Stop Solution	10 mL	1 bottle	Ready-to-use, 100 μ L/well
		ELISA Plate	/	1 plate	High affinity, removable plate
		Plate Sealers	/	4 strips	Ready-to-use
		Instruction	/	1 copy	Ready-to-use

* All reagents stored at 2 $^{\circ}$ C~8 $^{\circ}$ C, shelf life of 12 months.

[Instructions for use]

1. The product can meet the use of 1 ELISA plate.
2. Ready-to-use reagents are working solutions and can be used directly.
3. Dilute the concentrated solution into working solution according to the instructions before use.
4. The usage of each hole is based on actual demand, and the recommended usage is for reference only.
5. TMB prevents contamination, toxicity, and skin contact.
6. The termination solution contains sulfuric acid and is corrosive. Avoid skin contact.
7. A detection wavelength of 450 nm and a reference wavelength of 610 nm~630 nm are recommended for post-termination readings.

[Company Information]

Manufacturer and after-sales service unit Name: Shenzhen Dakewe Bio-engineering Co., Ltd.

Website: www.dakewe.com

Telephone: (86-755) 86235300

Email: RD@dakewe.com

Address: Room 702-703, Building No.1, Shenzhen Biomedicine Innovations Industrial Park, No.14 Jinhui Road, Kengzi Street, Pingshan District, Shenzhen, China

After-sales service telephone: (86-755) 86235300

Zip Code: 518122